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Large Capacity Refrigerated Centrifuge
for blood bags

9942

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Large Capacity Refrigerated Centrifuge

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Large Capacity Refrigerated Centrifuge for blood bags

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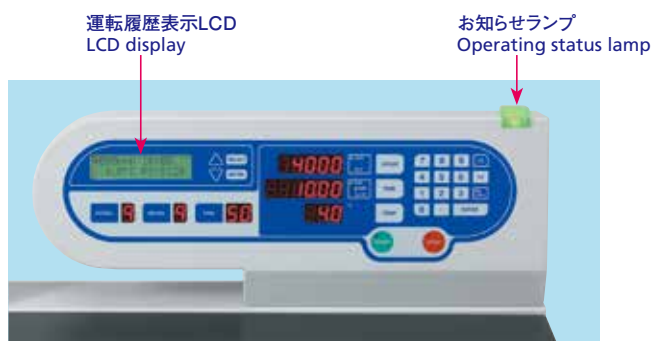
Counterbalanced lid opens and closes easily, with safety provided by the electrical interlock system.
A single Model 9942 unit can centrifuge 12 blood bags at once.

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LCD display functions:

- Setting operation parameters.
Number of cycles of use/Time of use/Error history.
Lamp color indicating stopped rotor.
Clock function (for LOG printer)
- Rotor parameter settings.
Max speed/RCF indications and changing of rotor radius.
Number of cycles of rotor use and rotor use time indications.

- An electrical interlock system is employed.
The chamber lid locks automatically and can be closed with a gentle push.
- Operating status lamp and buzzer indications let you know what the centrifuge is doing at a distance.
 - Seven different lamp colors can be selected as desired for each program number, to indicate when the centrifuge is running or stopped.
 - Five unique buzzer melodies, beeping, or silent mode can be selected.
- The inner cup for blood bags is designed for ease of use.
Special bottom surface design facilitates removing blood bags with minimal tilting of the inner cup, to preserve centrifuged layering.
- Acceptable imbalance up to 50 grams.
Measured and collected blood bags can be centrifuged directly. If imbalance is excessive, an imbalance sensor automatically stops the centrifuge.
- Maximal reproducibility of separation states.
Integrator g-sec values (integrated centrifugal values) can be measured and set.
- Step centrifugal function is available.
This centrifuge is compatible with automatic umbilical cord blood isolation unit.
- Complies with GMP requirements.
 - Easy to check actual rotor speed with the tachometer port.
 - LOG system and bar-code reader are available (optional).

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⚠️ ロータRS-8120には耐用年数と耐用回数があります。取扱説明書をご覧ください。
 Rotor RS-8120 have prescribed usage lifetimes (durable period) and durable cycles, after which they must be replaced. Please refer to the Instruction & Service Manual.

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- 400/500mL blood bags are inserted in the inner cup that has a central separator.
The separator prevents bag contamination if a bag breaks during operation.
- This inner cup can be used with both RS-8120 (for model 9942) and RS-8080 (for model 9920) rotors.
(Code No.055-6226/6pcs/set and No.055-6246/6pcs/set for RS-8120, Code No.055-6220/4pcs/set and No.055-6244/4pcs/set for RS-8080).
- Adaptor for 200mL blood bags is also available. (Code No.055-6290/12pcs/set for RS-8120, Code No.055-6298/8pcs/set for RS-8080).
- Bottom surface is designed to minimise tilting when removing blood bags from one side.

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- In the use of certain commercially available inner cups, separated layers are often disturbed when removing bags in one side while bags in the other side kept inside.
In such cases, the inner cup must be placed on a special board and disturbed blood bags must be re-centrifuged, requiring additional time and cost.

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Large Capacity Refrigerated Centrifuge

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Easily viewable LED display and LCD display for setting/error indication.
6 x 1L large volume high-speed angled rotor.

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*1 1,000mL Herolab bottles (nominal capacity) (Catalogue No.253562, 253565/polycarbonate, No.253572, No.253575/polypropylene), 1,000mL Nalgene bottles (nominal capacity) (Catalogue No.3120-1000/polypropylene copolymer, No.3122-1000/polycarbonate) are available.

*2 Sealing caps are equipped with bottle (Code No.052-1530/2pcs).

*3 Sealing caps are equipped with bottle (Code No.052-7080/2pcs).

*4 1,000mL Nalgene bottles (actual capacity) (Catalogue No.3140-1006/polycarbonate, No.3141-1006/polypropylene) are available.

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⚠️ Rotors RS-6602, RS-7002, RS-7100 and RS-8080 have prescribed usage lifetimes (durable period) and durable cycles, after which they must be replaced.
 Please refer to the Instruction & Service Manual.

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*1 This tube rack is a two-compartment system and easy to carry.
 *2 Equipped with a custom ordered adaptor. The adaptor is a custom order (charged option). Please let us know the type of rack you intend to use. Other adaptors can be made. Please contact your local distributor.
 *3 Custom ordered tube racks can be made. Please contact your local distributor.
 *4 Please contact your local distributor before ordering K12510 fixed-angle rotor. Sealing caps are equipped with 1,000mL stainless bottle. With rotor lid, 6,000rpm or less (9920), 6600rpm (9940), with sample temperature below 4 degrees.
 *5 Polycarbonate and polypropylene bottles are available.

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⚠️ RS-3900, RS-3920 and K12510 rotors have prescribed usage lifetimes (durable period) and durable cycles, after which they must be replaced. Please refer to the Instruction & Service Manual.

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(*) 1) Not including projections.

* Use only rotors, adaptors and tube racks made by KUBOTA. Do not attempt to use rotors, adaptors or tube racks made by other manufacturers.

* Follow all instructions in the relevant Instruction & Service Manual when operating centrifuges.

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URL <http://www.centrifuge.jp>

Kubota has acquired ISO 9001 and ISO 13485 certification.

NOTE

Products in this catalogue are designed for use by people having the requisite technical knowledge, and must always be used with care. People who do not have adequate technical knowledge or training should use our products only under appropriate supervision, or else hazards are likely to occur.

DISCONTINUE USE INDICATIONS

Always discontinue use if any of the conditions listed at right are evident.

- The rotor or bucket appears to be damaged, cracked or corroded.
- The rotor reaches the time for replacement (durable years) or durable cycles.
- The equipment emits a burnt smell.
- You receive a minor electric shock when you touch the equipment with your bare hand.
- Any other abnormality or indication of failure is noticed.

IMPORTANT

In the event of any observed abnormality in the equipment or its parts, to prevent occurrence of functional failure or accidents, immediately turn off the power switch, disconnect the power cable plug from the mains socket and contact your local distributor or Kubota Corporation if there is no local distributor in your neighbourhood.

Safety Precautions

For your safe use of the equipment, be sure to read the instruction manual carefully before you start operations.

Do not misplace the instruction manual. Keep the instruction manual nearby so you can refer to it whenever necessary.

- The term of supplying spare parts for repair is 10 years after discontinuation of production (except spare parts which we are unable to procure).
- This catalogue is not for distribution in the USA, Canada and Mexico as products shown are not for sale in these countries.

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