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Large Capacity Refrigerated Centrifuge
for blood bags

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Large Capacity Refrigerated Centrifuge for blood bags

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Counterbalanced lid opens and closes easily, with safety provided by the electrical interlock system.
A single Model 9942 unit can centrifuge 12 blood bags at once.

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LCD display functions:

- Setting operation parameters.
Number of cycles of use/Time of use/Error history.
Lamp color indicating stopped rotor.
Clock function (for LOG printer)
- Rotor parameter settings.
Max speed/RCF indications and changing of rotor radius.
Number of cycles of rotor use and rotor use time indications.

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■ An electrical interlock system is employed.

The chamber lid locks automatically and can be closed with a gentle push.

■ Operating status lamp and buzzer indications let you know what the centrifuge is doing at a distance.

- Seven different lamp colors can be selected as desired for each program number, to indicate when the centrifuge is running or stopped.
- Five unique buzzer melodies, beeping, or silent mode can be selected.

■ The inner cup for blood bags is designed for ease of use.

Special bottom surface design facilitates removing blood bags with minimal tilting of the inner cup, to preserve centrifuged layering.

■ Acceptable imbalance up to 50 grams.

Measured and collected blood bags can be centrifuged directly. If imbalance is excessive, an imbalance sensor automatically stops the centrifuge.

■ Maximal reproducibility of separation states.

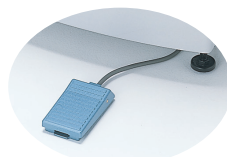
Integrator g·sec values (integrated centrifugal values) can be measured and set.

■ Step centrifugal function is available.

This centrifuge is compatible with automatic umbilical cord blood isolation unit.

■ Complies with GMP requirements.

- Easy to check actual rotor speed with the tachometer port.
- LOG system and bar-code reader are available (optional).

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RS-8120 Rotor especially designed for blood bags (only for model 9942)

Maximum Speed 4,000rpm
 Maximum RCF 5,240xg
 Maximum Capacity 血液バッグ x 12 12 x Blood bag

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Rotor RS-8120 have prescribed operating lifetime and years of operation, after which they must be replaced.
 Please contact us for more information.

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- In the use of certain commercially available inner cups, separated layers are often disturbed when removing bags in one side while bags in the other side kept inside.
 In such cases, the inner cup must be placed on a special board and disturbed blood bags must be re-centrifuged, requiring additional time and cost.

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Large Capacity Refrigerated Centrifuge

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Easily viewable LED display and LCD display for setting/error indication.
6 x 1,000mL large volume high-speed angled rotor.

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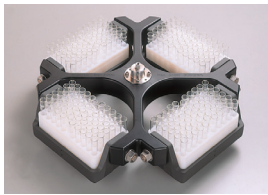
⚠ Rotors RS-6602, RS-7002, RS-7100 and RS-8080 have prescribed operating lifetime and years of operation, after which they must be replaced. Please contact us for more information.

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
*1 1,000mL Herolab bottles (Catalogue No.253562, 253565/polypropylene, No.253572, 253575/ polycarbonate), 1,000mL Nalgene bottles (Catalogue No.3120-1000/polypropylene copolymer, No.3122-1000, No.3140-1006/polycarbonate, No.3141-1006/polypropylene) are available.
*2 Stainless bottle (code No.052-1530/2 pcs) are available.
*3 Stainless bottle (code No.052-7080/2 pcs) are available.
*4 Code No.055-6220 is adapter for inner cup (code No.055-6220).

9920/9940 ロータ表 Rotor Table for the model 9920/9940

RS-3900 (only for model 9920)
Maximum Speed 3,000rpm
Maximum RCF 2,490xg
Maximum Capacity 3,780mL

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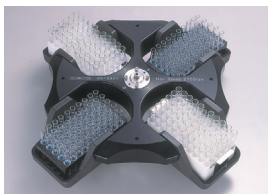RS-3920 (only for model 9940)
Maximum Speed 2,700rpm
Maximum RCF 2,140xg
Maximum Capacity 3,900mL

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⚠️ RS-3900, RS-3920 and K12510 rotors have prescribed operating lifetime and years of operation, after which they must be replaced. Please contact us for more information.

- *1 The adaptor is a custom order (charged option). Please let us know the type of rack you intend to use.
- *2 Custom ordered tube racks can be made. Please contact your local distributor.
- *3 Stainless bottle (code No.K13510/6 pcs) are available.
- *4 1,000mL Herolab bottles (Catalogue No.253562, 253565/polypropylene, No.253572, 253575/polycarbonate) are available.

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Kubota has acquired ISO 9001 and ISO 13485 certification.

Precautions for use Products in this catalogue are designed for use only by people who have the requisite technical knowledge, and must always be used with considerable care and only for their intended purpose. People who do not have adequate technical knowledge or training should only use the products under appropriate supervision by someone with expert knowledge, or else accidents are likely to occur.

Prohibited Please immediately stop using the products in any of the cases listed on the right.

- The rotor or buckets appear to be damaged or corroded.
- When the replacement period (years of operation, operating lifetime) of a rotor has passed.
- The equipment emits a burning smell or becomes abnormally hot.
- You receive a weak electric shock when you touch the equipment with your bare hands.
- When any other abnormality or indication of failure is noticed.

IMPORTANT If any of the cases listed at the left occurs, immediately turn off the power, disconnect the power cable plug or connecting terminals from the main power outlet, place a "Do not use" sign on the unit, and contact the nearest branch of Kubota Corporation.

Safety Precautions To use the equipment safely, be sure to read the instruction manual carefully before you start operations.

Do not misplace the instruction manual. Keep the instruction manual nearby so that you can refer to it whenever necessary.

- The term of supplying spare parts for repair is 7 years after discontinuation of production (except spare parts which we are unable to procure).
- This catalogue is not for distribution in the USA, Canada and Mexico as products shown are not for sale in these countries.

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